

Hydros Unpacking Instructions

Step 01

Remove the packing tape and open the box.



Step 02

Remove tray of accessories from the top of the box.



Step 03

Put box vertically and slide box off of the Hydros deioniser.



Step 04

Remove Styrofoam and plastic bag from the Hydros Deioniser.



Step 05

Hydros deioniser is now unboxed.





Contact Us

For more information, contact
Fistreem directly at:

Beacon House,
Nuffield Road,
Cambridge,
CB4 1TF, UK



+44 (0)1223 727 100



+44 (0)1223 727 101



support@fistreem.co.uk



www.fistreeminternational.com

