

How to replace your pretreatment cartridge

Step 01

Ensure the Inlet flow is turned off and that the remaining water has been drained out.



Step 02

Remove the two hose pipes connected to the cannister, found at the top and bottom.



Step 03

Put the old cartridge aside and place the new cartridge on the holder.



Step 04

Set up the layout of the cannister as shown, with the white valve inserted on the cannister and the grey connector at the bottom. You will have to unscrew the end caps slightly on the valve and connector to allow for it to slide into the nozzle.



Step 05

Screw the end caps of the valve and connector to make a tight seal.



**Step
06**

Remove the other end cap of the connector.

**Step
07**

Attach the inlet hose (the hose connected to the tap) to the connector.

**Step
08**

Remove the top end cap of the valve.

**Step
09**

Attach the reinforced hose pipe to the valve.

**Step
10**

Attach the other end of the reinforced pipe into the Pre-Treated Water socket located at the left side of the Still marked with the label shown in the image or to where your system connects to.

