

ProcScan - Setting Up Quick Guide

1 Install scanner software

- ➔ Install the scanner software onto ProtoCOL 3 or PC according to scanner instructions
- ➔ Connect USB cable to ProtoCOL 3 or PC

2 Start ProtoCOL 3 software



- ➔ Log on to ProtoCOL 3 software

3 Twain source


Click



Select



Select

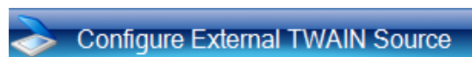

- ➔ Click on the ProtoCOL 3 icon in the top left of the screen
- ➔ From the drop down menu select Settings then "Configure External TWAIN Source"
- ➔ Select the scanner being used

4 Select scanner

Click



Click



- ➔ Select the Image tab
- ➔ Select the scanner by selecting TWAIN Device from the drop down menu on the left side of the screen labelled Internal Camera if scanner is being used with a ProtoCOL 3

5 Capture image

Click



- ➔ Click Capture Image
- ➔ The scanner software will now open and the settings can be set for the type of plate to be analysed

CONTACT SYNBIOSIS:
EUROPE:
Tel: +44 (0)1223 727125
Email: support@synbiosis.com

USA:
Tel: 800 686 4451/301 662 2863
Email: support@synbiosis.com

www.synbiosis.com
PSSUQ60212