

Protos 3 Software Update Quick Guide

1 Start PC

- ➔ Log onto PC as an Administrator and shut all running software down


2 Insert software source




- ➔ Insert the software source, unzipping if the installation set is zipped. Run the file Setup.exe and follow the on-screen instructions.

The PC may require a Windows update (particularly if the PC is Windows XP, Windows Vista or Windows 7). If you are presented with a message stating "**This application needs Microsoft .Net Framework 4.0 full version which is not yet installed. This can be found in the Support directory.**", then follow step 3 if you have a full installation set or 4 if you just have an update set.

3 Support folder

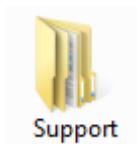

- ➔ Open the Support folder

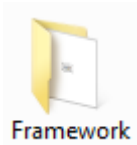

- ➔ Open the Framework folder



- ➔ Run SetupFramework.exe. Follow the on-screen instructions, performing a system restart if update required.



- ➔ Run Setup.exe once again and follow the on-screen instructions (see Protos 3 Full Installation Guide for details).

4

Download



dotNetFx40_Full_x86_x64.exe



Open a web browser, download and run the .NET Framework 4.0 full update from:

<http://www.microsoft.com/en-gb/download/details.aspx?id=17718>

5



Setup.exe



Run Setup.exe once again and follow the on-screen instructions (see Protos 3 Full Installation Guide for details).

Protos 3 is now ready to be used

CONTACT SYNBIOSIS:
EUROPE:
Tel: +44 (0)1223 727125
Email: support@synbiosis.com

USA:
Tel: 800 686 4451/301 662 2863
Email: support@synbiosis.com

www.synbiosis.com
P3si0614