

Condenser Foam Implementation Guide for Fistreem III 4L/Hr Still


Step 01

Ensure the still is turned off and has cooled down then remove the front panel.



Step 02

Find the foam in the accessory pack and peel the paper off the tape carefully (don't tear the foam) and slightly squash the foam to fit.

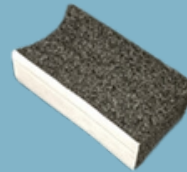

Step 03

Slide the foam up into the top mold and in front of the condenser whilst holding on to the condenser to avoid the glass from moving.



Step 04

Stick the foam on the inside of the top mold ensuring it is aligned to the face of the condenser.



Step 05

Replace the front panel.

