



ProcScan Unit set-Up & Modification

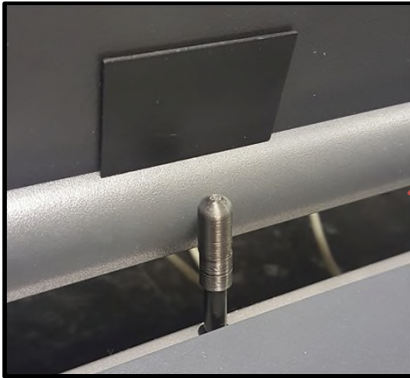



Place the black cover over the square hold in the CPU and place the extender piece over the peg to ensure it is adequately pushed down. If the peg is not pushed down the scanner will not work

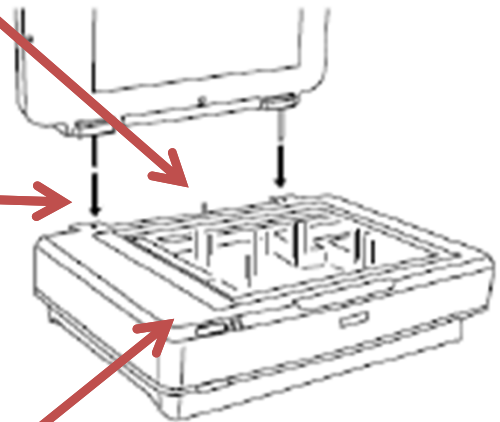

Place the spacer over the prongs for the TPU on the scanner lid




Stick the spacers on the front of the scanner body as shown

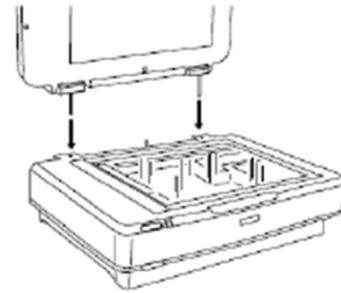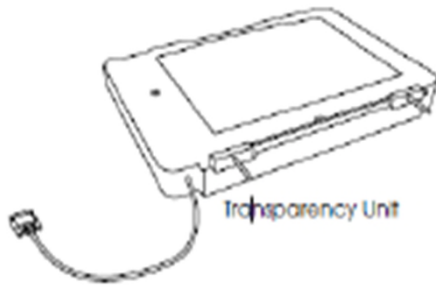

Attach Transparency unit to the top of the scanner and plug in connection wires at the back of the unit

Unlock the scanner arm within the Transparency unit by twisting it to the Unlock position

Use a coin to unlock the transportation lock by turning it counterclockwise one-quarter.

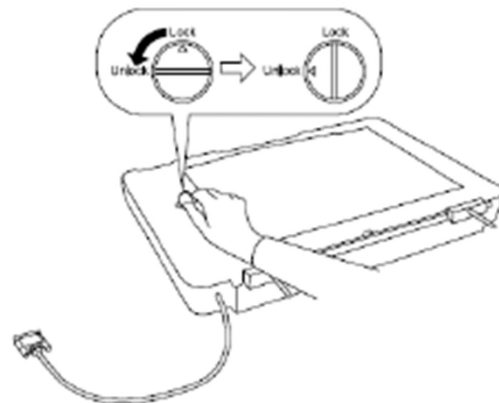

Note: if this is not done the Transparency unit will not function

Your ProcScan is now ready to use with your PC or ProtoCOL 3 system.

CONTACT SYNBIOSIS:

EUROPE:
Tel: +44 (0)1223 727125
2863
Email: support@synbiosis.com

USA:
Tel: 800 686 4451/301 662
Email: support@synbiosis.com

www.synbiosis.com