

# ProtoCOL 3 Software Update Quick Guide

## 1 Start PC

- ➔ Log onto PC as an Administrator and shut all running software down


## 2 Insert software source




- ➔ Insert the software source, unzipping if the installation set is zipped. Run the file Setup.exe and follow the on-screen instructions.

The PC may require a Windows update (particularly if the PC is Windows XP, Windows Vista or Windows 7). If you are presented with a message stating "**This application needs Microsoft .Net Framework 4.0 full version which is not yet installed. This can be found in the Support directory.**", then follow step 3 if you have a full installation set or 4 if you just have an update set.

## 3 Support folder

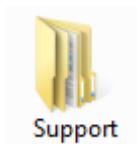

- ➔ Open the Support folder

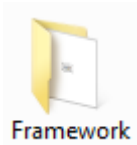

- ➔ Open the Framework folder



- ➔ Run SetupFramework.exe. Follow the on-screen instructions, performing a system restart if update required.



- ➔ Run Setup.exe once again and follow the on-screen instructions (see ProtoCOL 3 Full Installation Guide for details within Installation -> Manuals-> Quick Guides).

4

## Download



dotNetFx40\_Full\_x86\_x64.exe



Open a web browser, download and run the .NET Framework 4.0 full update from:

<http://www.microsoft.com/en-gb/download/details.aspx?id=17718>

5



Setup.exe



Run Setup.exe once again and follow the on-screen instructions (see ProtoCOL 3 Full Installation Guide for details within Installation -> Manuals-> Quick Guides).

**ProtoCOL 3 is now ready to be used**

CONTACT SYNBIOSIS:  
EUROPE:  
Tel: +44 (0)1223 727125  
Email: [support@synbiosis.com](mailto:support@synbiosis.com)

USA:  
Tel: 800 686 4451/301 662 2863  
Email: [support@synbiosis.com](mailto:support@synbiosis.com)

[www.synbiosis.com](http://www.synbiosis.com)  
P3si0614