

# ProtoCOL 3 Plus

## Connecting Touch Screen PC - Quick Guide


1

### Unpack ProtoCOL 3

Unpack the instrument and the PC and place on a secure level surface

### Attach bracket and PC to ProtoCOL 3 body

2

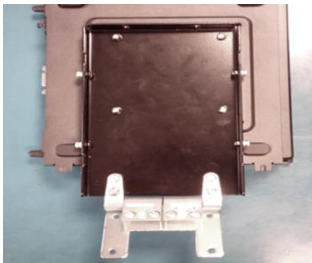

Attach the black bracket with PC to the instrument via the silver mounting plate as shown. The four screws are secured using a Pozidriv screwdriver (not supplied)



There are three leads that come out of the front of the ProtoCOL 3 body

### Attach monitor leads

3



Attach the 2 monitor leads to the monitor before attaching the monitor to the PC/bracket. Ensure that the ferrite (red arrow) is further from the monitor so that it does not impede monitor positioning

## Attach leads to PC and monitor

4



There are 3 leads that come from the front of the ProtoCOL 3 system. These will be labelled to explain where to connect each lead

Attach the leads to the PC and to the monitor before securing the monitor onto the bracket (step 6)

## Attach monitor to PC/bracket

5



Slide the bracket that is attached to the monitor over the top, lining the two pins with holes in monitor bracket - taking care to feed the cables under the monitor so as not to impede the movement of the monitor

Secure the monitor onto the bracket using the 4 screws

## Attach power cords to the mains power blocks

6



There are 2 power leads found at the back of the ProtoCOL 3 system. Attach the mains cables provided to the power blocks and plug into a socket

THE LOGIN DETAILS FOR THE PROTOCOL 3 PLUS  
ARE

**Username: P3Admin**

**Password: The unit's serial number in capitals**



**Your ProtoCOL 3 Plus is now ready to use!**

### CONTACT SYNBIOSIS:

EUROPE:

Tel: +44 (0)1223 727125

Email: [support@synbiosis.com](mailto:support@synbiosis.com)

[www.synbiosis.com](http://www.synbiosis.com)

USA:

Tel: 800 686 4451/301 662 2863

Email: [support@synbiosis.com](mailto:support@synbiosis.com)

P3CTSPC0618